

What's New?

Software Update | July 31, 2018

1. Languages

- a. Added new languages: German, Chinese (Simple), Chinese (Traditional), Spanish, and Italian along with the existing French (Canadian), French (European), and English.
- b. When a language other than English is selected, and the Quantifit is connected but has not been calibrated, the user will have an option to view calibration instructions in the selected language.

2. Installation

- a. Installation no longer requires a re-boot.
- b. Dot-Net Framework set automatically during Windows 10 install.

3. Import/Export

- a. Updated personnel import template use Excel .xlsx.
- b. A blank Personnel Import template is available for download on the Personnel Import screen.
- c. Automatic file correction when importing data from the Quantifit.

4. Settings

- a. Ability to change fit test expiration date to 2 years to comply with Canadian regulations. This 2-year expiration is visible when using a card printer.
- b. Predetermined fields selected by default on Company Edit screen.
- c. Ability to hide unused Protocols.

5. General Improvements

- a. Start Button on Backup function removed.
- b. Employee count displays at top of Personnel Screen (also works when filtering the employee list.)
- c. Backup reminder defaults to 0 days, which will ask operator to back up database each time the software is closed.
- d. Test screen will remember last protocol used.
- e. Software version viewable on the software title bar.
- f. If the Quantifit is attached after the software is started, the software will import the current calibration data. (It is recommended that a calibration is completed and Quantifit connected to computer before launching software.)